Free and Total Chlorine Test Kit

0–0.7 and 0–3.5 mg/L

To ensure accurate results, read carefully before proceeding.

WARNING: Handling chemical samples, standards, and reagents can be dangerous. Review the Material Safety Data Sheets before handling any chemicals.

General Test Information

- Wash all labware between tests. Contamination may alter test results. Clean with a non-abrasive detergent or a solvent such as isopropyl alcohol. Wipe and dry with a soft cloth. Do not use paper towels or tissue on plastic tubes as this may scratch them. Rinse with deionized water.
- High concentrations of monochloramine will interfere with Free Chlorine test results.

Using PermaChem® Powder Pillows:


1. Tap the bottom of the pillow on a hard surface.
2. Tear open the pillow along the dashed line.
3. Push the side edges inward to form a spout.
4. Pour the contents into the sample.

Measuring Hints

- Accuracy is not affected by undissolved powder.
- Rinse all viewing tubes thoroughly with sample before testing.
- For optimum test results, check reagent accuracy with each new lot of reagents. Use the standard solution listed in *OPTIONAL REAGENTS AND EQUIPMENT* below. Follow the instructions included with each standard solution.

- If the color from the Low-Range test reads off the high end of the scale, repeat the test using a new sample and the High-Range instructions.

Figure 1 Using the Long Path Viewing Adapter in the Color Comparator

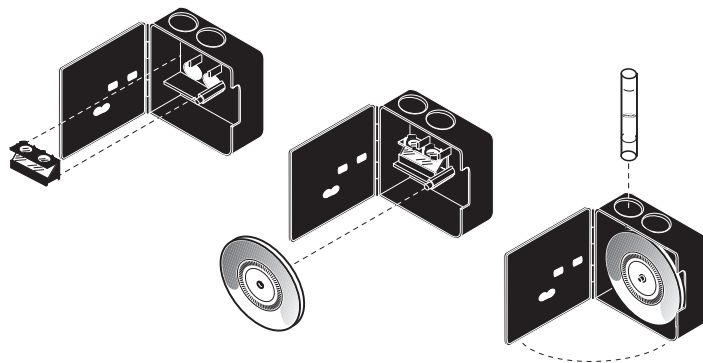Free Chlorine or Total Chlorine (Low Range) Test, 0–0.7 mg/L

1. Insert the Long Path Viewing Adapter into the color comparator (see Figure 1 above).
2. Fill a viewing tube to the top line with sample. This is the sample blank.
3. Place this tube in the top left opening of the color comparator.
4. Fill the square bottle to the 25-mL mark with sample.
5. Add the contents of one DPD Free Chlorine Reagent Powder Pillow or one DPD Total Chlorine Reagent Powder Pillow to the bottle. Swirl to mix.

Note: If testing Free Chlorine, complete the test and read the result within one minute of adding the reagent.

- Note:** If testing Total Chlorine, read the test result after three minutes but before six minutes of adding the reagent.
6. Fill another viewing tube to the top (15 mL) line with the prepared sample.
 7. Place this tube in the top right opening of the color comparator.
 8. Hold the comparator with the tube tops pointing toward a bright light and look through the openings in front.
 9. Rotate the color disc until the colors match.
 10. Divide the value in the scale window by 5 to obtain the mg/L free or total chlorine.

Free Chlorine or Total Chlorine (Mid Range) Test, 0–3.5 mg/L

1. If the color comparator has the Long Path Viewing Adapter in place, remove it (see Figure 1 above).
2. Fill a viewing tube to the bottom (5-mL) line with sample. This is the sample blank.
3. Place this tube in the top left opening of the color comparator.
4. Fill the square bottle to the 25-mL mark with sample.
5. Add the contents of one DPD Free Chlorine Reagent Powder Pillow or one DPD Total Chlorine Reagent Powder Pillow to the bottle. Swirl to mix.

Note: If testing Free Chlorine, complete the test and read the result within one minute of adding the reagent.

- Note:** If testing Total Chlorine, read the test result after three minutes but before six minutes of adding the reagent.
6. Fill another viewing tube to the bottom (5-mL) line with the prepared sample.
 7. Place this tube in the top right opening of the color comparator.
 8. Hold comparator up to a bright light and look through the openings in front.
 9. Rotate the color disc until the colors match.
 10. Read the mg/L free or total chlorine in the scale window.

REPLACEMENT PARTS

Description	Unit	Cat. No.
Bottle, square w/ 25-mL mark.....	each	23510-00
Color Comparator Box	each	1732-00
Color Disc, DPD Chlorine, 0–3.5 mg/L	each	21988-00
Color Viewing Tube, plastic, with cap	4/pkg	46600-04
DPD Free Chlorine Reagent Powder Pillows.....	100/pkg	14070-99
DPD Total Chlorine Reagent Powder Pillows.....	100/pkg	14064-99
Instruction Card, CN-70 Test Kit	each	14542-38
Long Path Viewing Adapter	each	24122-00

OPTIONAL REAGENTS AND EQUIPMENT

Description	Unit	Cat. No.
Caps, for Color Viewing Tubes 46600-04	4/pkg	46600-14
Chlorine Standard Solution, 50–75 mg/L, 2-mL PourRite™ Ampule	20/pkg	14268-20

PourRite™ and PermaChem® are trademarks or registered trademarks of Hach Company.

FOR TECHNICAL ASSISTANCE, PRICE INFORMATION AND ORDERING:

In the U.S.A. - Call toll-free 800-227-4224
 Outside the U.S.A. - Contact the HACH office or distributor serving you.
 On the Worldwide Web - www.hach.com; E-mail - techhelp@hach.com

HACH COMPANY
 WORLD HEADQUARTERS
 Telephone: (970) 669-3050
 FAX: (970) 669-2932

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



FOR TECHNICAL ASSISTANCE, PRICE INFORMATION AND ORDERING:
In the U.S.A. - Call toll-free 800-227-4224
Outside the U.S.A. - Contact the HACH office or distributor serving you.
On the Worldwide Web - www.hach.com; E-mail - techhelp@hach.com

HACH COMPANY
WORLD HEADQUARTERS
Telephone: (970) 669-3050
FAX: (970) 669-2932

Free and Total Chlorine Test Kit 0–3.5 mg/L

WARNING: *Handling chemical samples, standards, and reagents can be dangerous. Review the Material Safety Data Sheets before handling any chemicals.*

Note: *Remove the Long Path Viewing Adapter from the Color Comparator.*

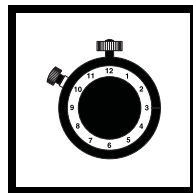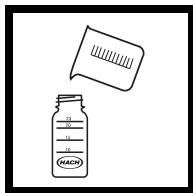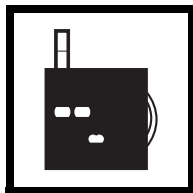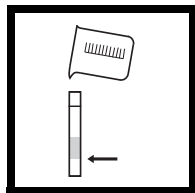English

1. Fill the tube to the bottom line with sample.

2. Place tube in the left opening of the color comparator.

3. Fill the square bottle to the 25-mL mark with sample.

4. Add contents of **appropriate** Powder Pillow to the bottle. Swirl.

Note: *Complete steps 5–9 within 1 min. of adding powder for Free Cl₂ and within 3–6 min. for Total Cl₂.*

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